

CALIBRATION CURVE

For accurate work, standard solutions should be included with each determination. Variations in chemical reactions of new batches of reagents and in instrument behavior combine to cause variability in the absorbance 'A' of samples. It is useful to document the successive absorbance 'A' readings of the standard for quality control purposes. Furthermore, when Beer-Lambert law is obeyed, calibration curve is prepared for the range of concentrations intended to be covered in practice. This curve is prepared by taking several dilutions of the substance to be measured, running the experiment as described, then plotting absorbance against concentration on a graph paper. When running the experiment, the concentration of the unknown is found from this curve.

Preparation of calibration curve

An alternative procedure to find the concentration of 'test' of unknown protein sample is to prepare a calibration curve, then read the concentration of the unknown.

Procedure

Three known concentrations of protein standard will be provided. Place 3 ml of each concentration in a glass tube and label (St₁, St₂ & St₃). Place 3 ml of distilled water in another tube and label B. Add 5 ml of Biuret colour reagent to all tubes, mix, incubate for 30 min. at room temp. or for 10 min. at 37 °C water bath. Read the absorbance of the standard and blank by the spectrophotometer at 540 nm against distilled water. Subtract 'B' reading from each reading of the standards, then construct a calibration or standard curve with absorbance on Y-axis and concentrations in g/L ml on X-axis as in fig. (1). Find the concentration of the unknown solution 'T' from the curve.

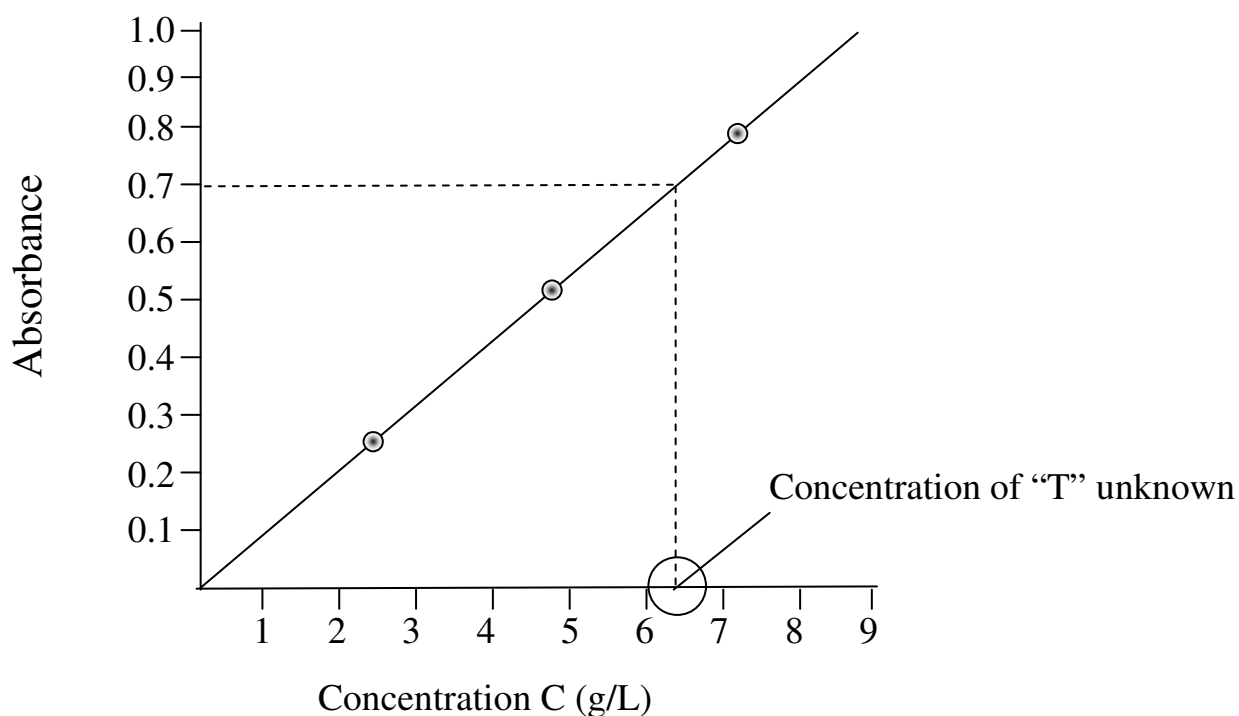


Fig. 1

$$C_T = \frac{A_T}{A_{St}} \times C_{St}$$

Reagents

- 1- Biurt reagent (as for total protein).
- 2- Protein standards : 2.5, 5.0, 7.5 g/L. They are equivalent to: 37.5, 75.0, 102.5 g/L in serum.

Prepare stock std: 5gm bovine albumin in 100 ml H₂O or (50 g/L).

- a. Take 12.5 ml of stock and dilute to 250 ml with DW. This will give 2.5 g/L.
 - b. Take 25 ml of stock and dilute to 250 ml with DW. This will give 5.0 g/L.
 - c. Take 37.5 ml of stock and dilute to 250 ml with DW. This will give 7.5 g/L.
- 3- Unknowns: the above standards can be given as unknowns or any other dilution could be made.